




## Maxi-Thermal Software

Works with c\temp-USB, edl-USB, 2c\temp-USB,  
and 2c\temp-RH-USB.



### Understanding the Menu Bar

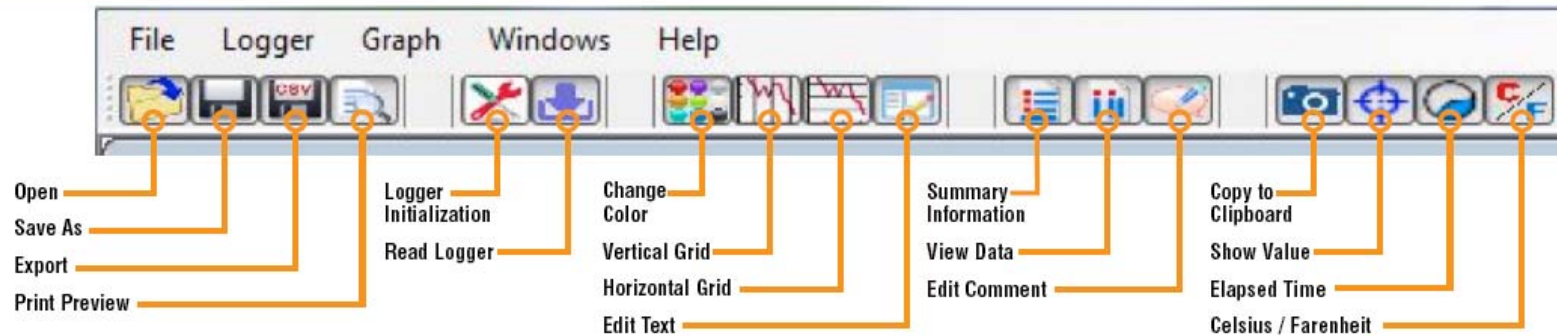

Q U I C K S T A R T Q U I D E

## Logger Initialization

1. Select Logger in menu bar
2. If there was a previous recording click **Erase All Data**
3. Enter **Description** for the recording
4. Select **Time Zone**
5. Select **Start Delay**
6. Select **Duration of Recording**
7. Select Units – Fahrenheit or Celsius
8. Select **No Alarm**, OR
9. Set **Maximum** (High Alarm) Temperature
10. Set **Minimum** (Low Alarm) Temperature
11. Click **Write to Device**

Logger Initialization

Description of Recording

Model number CTempUSB(Type E)

Description Maxitrak

(Max. 20 characters and can't have <>?/:\"/>

Current Clock

(GMT-08:00) Pacific Time (US & Canada)

3/6/2012 14:00:09

Logger Clock

(GMT-08:00) Pacific Time (US & Canada)

3/6/2012 14:00:09

Recording

Start Delay 5 0 Min.

Duration 6 60 Days

(Sample Interval: 10 Min 38 Sec)

Unit

☐ °F ☒ °C

☒ No alarm

Alarm

Maximum 30 °C

Minimum 20 °C

Quarantine 0

Event Log:

14:00:03 Found: cTemp\_usb on 682b4456d9&0#{53}

Write to device


Erase All Data

Extra settings...


SCREENSHOT 2 – Logger Initialization

## Read Logger


1. Select Read Logger on menu bar  
If the logger is recording you will see SCREENSHOT 3.
2. Select **Yes** or **No**
3. If you select Yes MaxiThermal will read the logger and display the graphical data (SCREENSHOT 4).

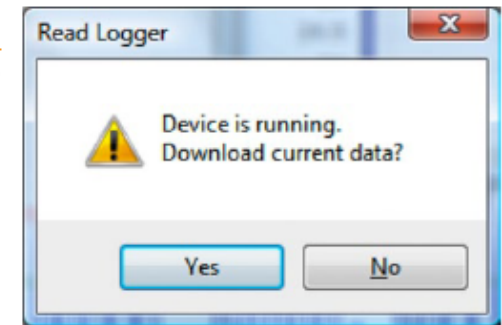


SCREENSHOT 3 – Read Logger

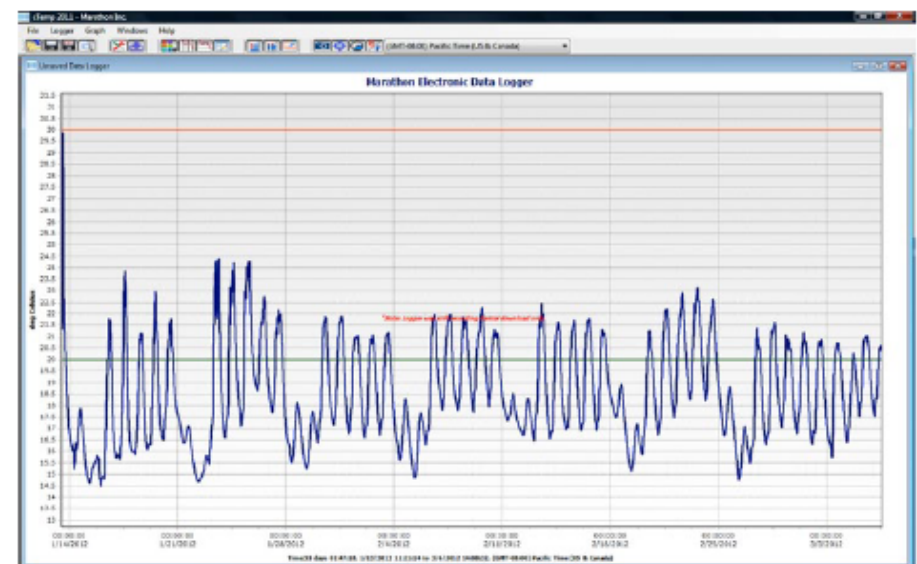


SCREENSHOT 4 – Graphical Data

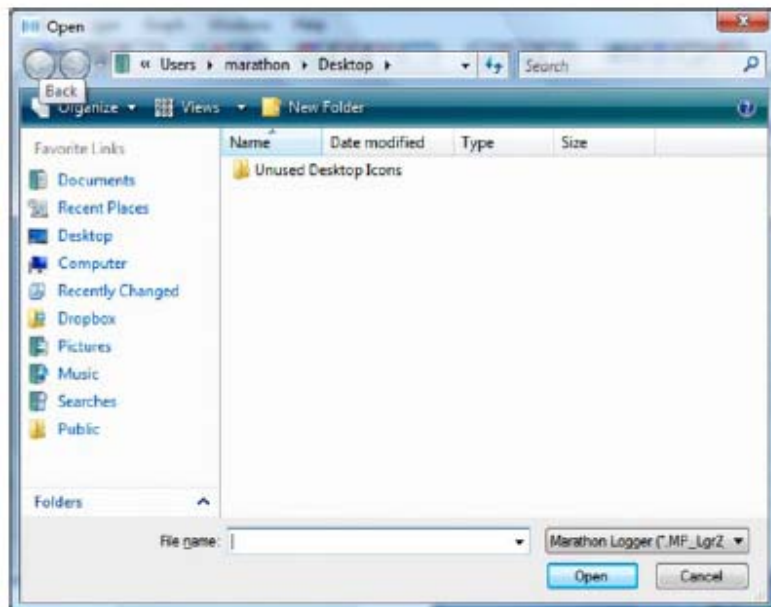


SCREENSHOT 6 – Open

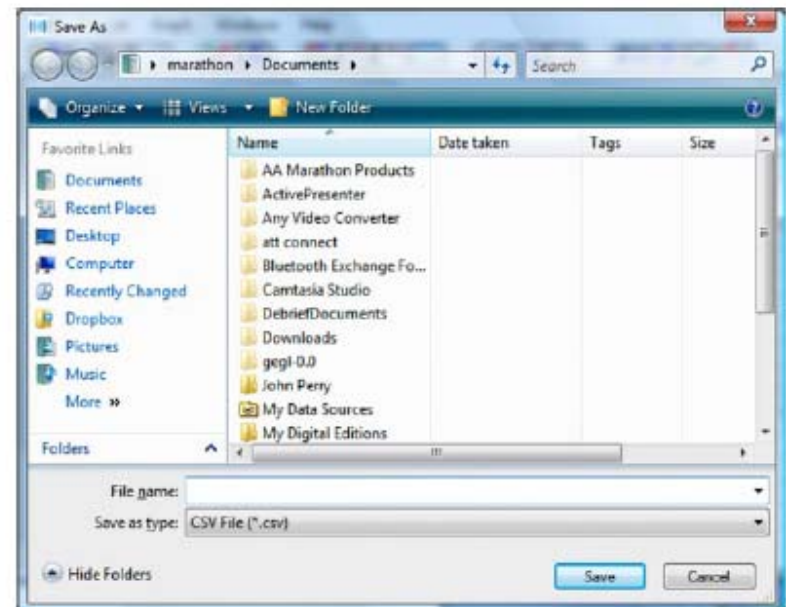


SCREENSHOT 8 – Export

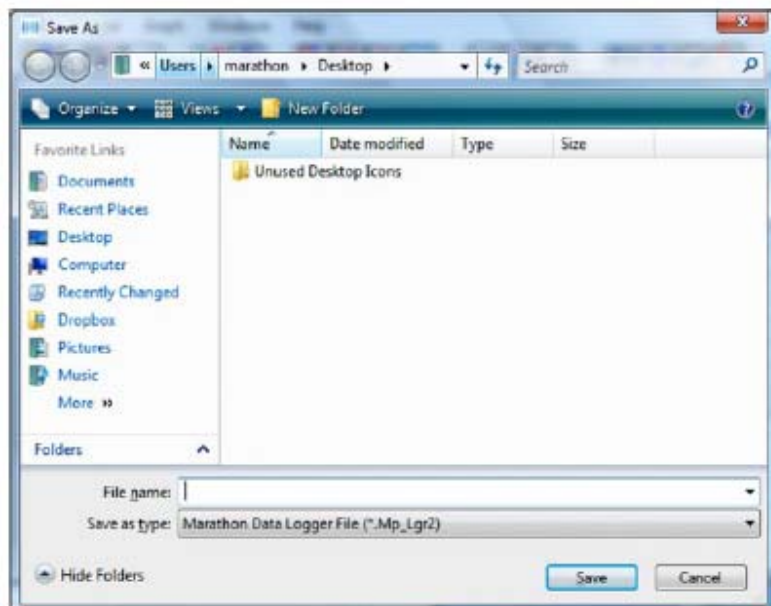

SCREENSHOT 7 – Save As

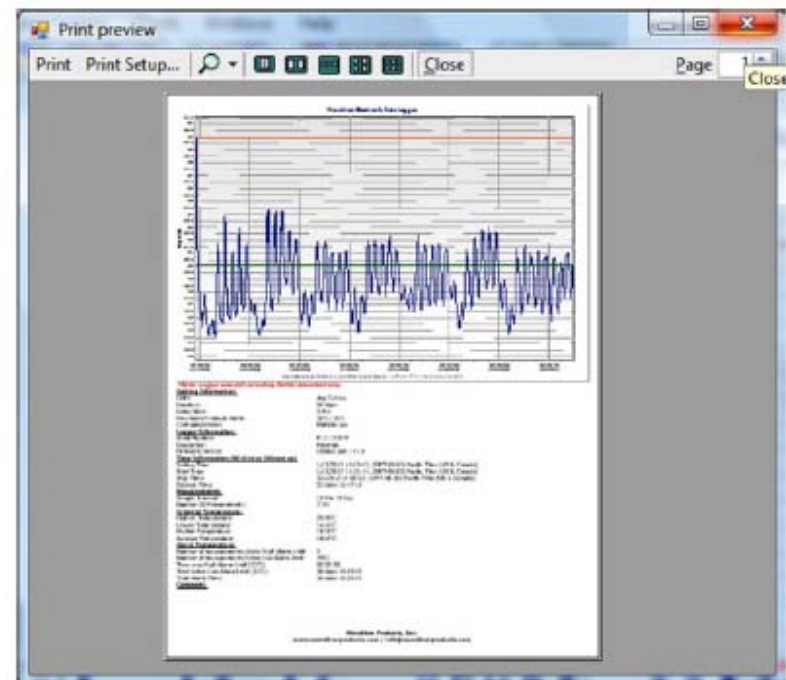

SCREENSHOT 9 – Print Preview

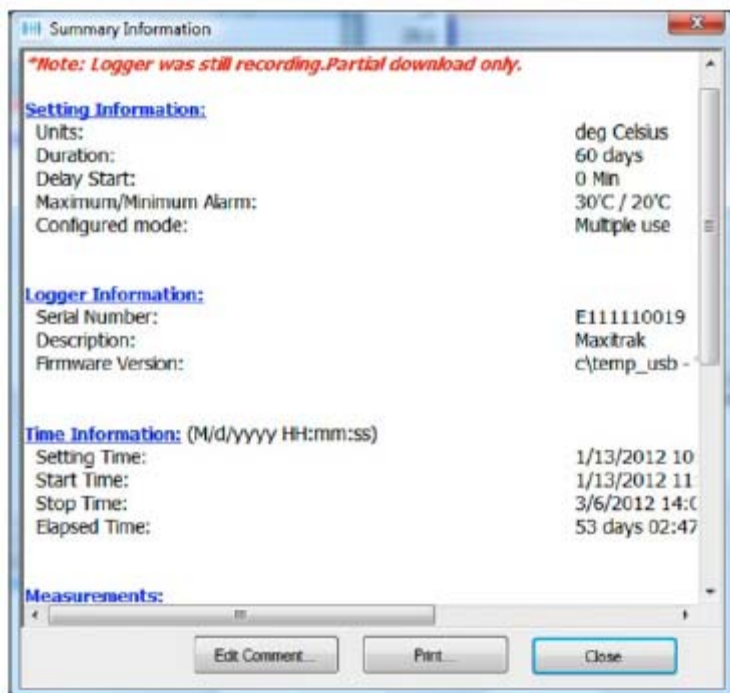

SCREENSHOT 10 – Summary Information

**View Data**

*\*Note: Logger was still recording. Partial download only.*

| TIME               | TEMPERATURE |
|--------------------|-------------|
| 1/13/2012 11:31:52 | 21.96 °C    |
| 1/13/2012 11:42:30 | 22.20 °C    |
| 1/13/2012 11:53:08 | 21.54 °C    |
| 1/13/2012 12:03:46 | 21.31 °C    |
| 1/13/2012 12:14:24 | 25.88 °C    |
| 1/13/2012 12:25:02 | 28.60 °C    |
| 1/13/2012 12:35:40 | 29.48 °C    |
| 1/13/2012 12:46:18 | 29.79 °C    |
| 1/13/2012 12:56:56 | 29.89 °C    |
| 1/13/2012 13:07:34 | 26.16 °C    |
| 1/13/2012 13:18:12 | 25.77 °C    |
| 1/13/2012 13:28:50 | 26.19 °C    |
| 1/13/2012 13:39:28 | 26.87 °C    |
| 1/13/2012 13:50:06 | 27.58 °C    |
| 1/13/2012 14:00:44 | 28.35 °C    |
| 1/13/2012 14:11:22 | 24.40 °C    |
| 1/13/2012 14:22:00 | 23.14 °C    |
| 1/13/2012 14:32:38 | 22.60 °C    |
| 1/13/2012 14:43:16 | 22.33 °C    |
| 1/13/2012 14:53:54 | 22.22 °C    |
| 1/13/2012 15:04:32 | 22.22 °C    |

Buttons: Print..., Close

SCREENSHOT 11 – View Data



1295 Morningside Ave Units 16, 17, & 18  
 Toronto ON M1B 4Z4 Canada  
 Telephone: 416-261-4865 Fax: 416-261-7879  
[www.scigiene.com](http://www.scigiene.com)