




	Value knob
	Function knob
	aspirating
	apocenososis
	Loud speaking symbol
	Power symbol



3. Operation

3.1 Power on



Keep pressing the value knob  2 seconds



3.2 Liquid handling

3.2.1 value selection



Left-right rotate the knob  quickly to active the volume adjusting mode, the “PI” symbol and volume value will be flashing under this mode. Rotate the knob  to do the value adjusting.

3.2.2 Aspiration & Apocenososis




 the arrow up symbol flashing when the dPette Electronic Pipettor is in aspirating mode, clicking the knob  to do aspirating.

 the arrow down symbol flashing when dPette Electronic Pipettor is in apocenososis mode, clicking the knob  to do apocenososis.




3.2.3 Mixing

Keep pressing the knob  under the aspiration mode will active the mixing mode. Under this mode, the dPette Electronic Pipettor will do the repetitive work of aspirating and apocenososis by the maximum speed 300uL. The maximum volume is 300-l in this function, If the setting volume is less than 300-l, the dPette Electronic Pipettor will do the mixing work by the set value. Release the knob , the mixture work will stop.




3.3 Function Setting

left-right rotate the function knob  quickly will active the function setting mode, rotate the function knob  to do function item selection; double click the value knob  to quit the function selecting mode.




3.1 Loud Speaker

Rotate the function knob  and select the loud speaker  symbol. Rotate the value knob  to do selection between “on” and “off”

3.2 Speed Adjusting

Rotate the function knob  until the arrow  symbol flashing. Rotate the value knob  to do the speed value adjusting.

3.3 Power Off



Rotate the function knob  to the power symbol  Press the value knob  to turn off dPette Electronic Pipettor.

dPette Electronic Pipettor (Part #: PPD01-24XX)



1295 Morningside Avenue, Unit 16-18
Scarborough, ON M1B 4Z4 Canada
Phone: 416-261-4865 Fax: 416-261-7879
www.scigiene.com

1. Structure

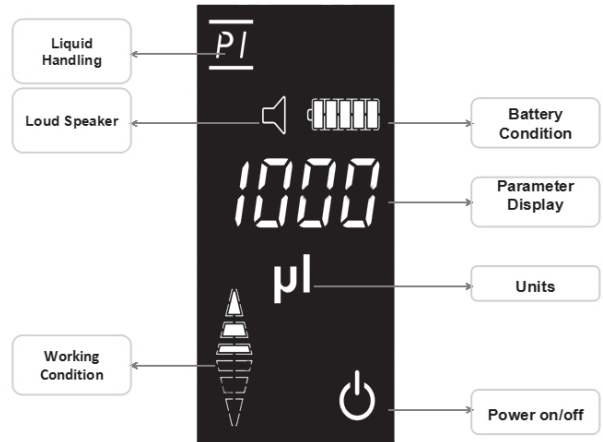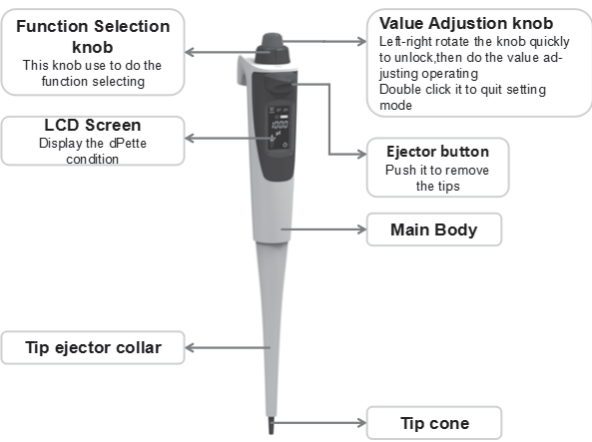

2. Charging and Reset

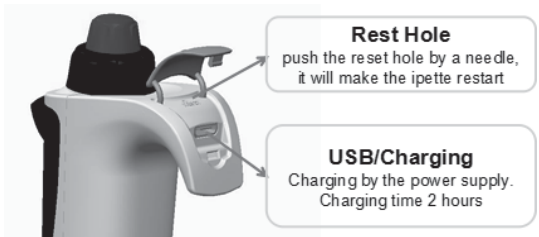


Image 3

